

OPERATING INSTRUCTIONS SOMNI CATCH



1. Position the SOMNI Catch under the stereotaxic nosecone.
2. Connect the active waste gas system (such as Somni EPS-3) tubing to the blue tubing adapter.
3. Connect the stereotaxic nosecone scavenging port to the black tubing already connected to the Catch.
 - a. Note: Choose correct tubing size based on manufacture of nose cone.
4. Verify the active waste gas system is turned on prior to using inhalant anaesthetic.
5. Attenuate the active waste gas scavenging system to at least 15 LPM negative flow.
6. At completion of inhalant anaesthesia, continue to use the active waste gas scavenging system for several seconds after the anaesthetic has been turned off to allow any remaining anaesthetic gas to be removed.

Cleaning: Remove the tubing attachments prior to cleaning with dilute chlorhexidine or similar disinfectant. Rinse and dry thoroughly.

SOMNI PROVIDES UNPARALLELED CUSTOMER SERVICE, CLINICAL AND TECHNICAL SUPPORT.

For more information or pricing, please contact our sales team.

☎ (T) 0800 052 5600 (D) 01872 248 890 (M) 07798969805

✉ enquiries@somniscientific.com 🌐 www.somniscientific.co.uk